

## The New MA160 Moisture Analyzer Manages your Sophisticated Tasks



# Fast, Precise and Easy Moisture Analysis

To be able to perform fast and precise analysis on a diverse range of product samples, you need a reliable instrument that gives you maximum flexibility.

The MA160 uses the thermogravimetric method to determine the moisture content of liquid, pasty and solid substances – conveniently, reliably and in a minimum of time. It delivers prompt, repeatable results and supports the development of new methods – in three uncomplicated and intuitive steps. The MA160 manages the generated methods professionally and securely and allows them to be transferred to other instruments.

Effortless cleaning

Fast and precise method development

High-speed measurements

Reliable performance testing

User-friendly operation

Large memory capacity

Display of measurement status



During a measurement, a status lamp indicates the current measurement status. Additionally, the ergonomic BetterClean design of the MA160 allows efficient and simple cleaning of the device.

Using integrated performance testing, the MA160 verifies its functionality at the touch of a button – ensuring permanent and flawless performance.





## Applications

With its optimized heating elements, the MA160 produces high-speed measurements. That's how the MA160's weighing system guarantees the required precision.


The MA160 is ideal for the moisture analysis of a varying range of samples under different conditions. It can generate new methods which allows you to create and effectively manage proprietary measurement procedures for various samples. This supports your work in the QC lab or in process control.

Typical areas of applications for the MA160 include the moisture analysis of foods, beverages, pharmaceuticals, chemicals, paper materials and environmental protection.

## Features

### High-speed measurements

Both the high-performance AURI heater and the sample chamber's geometry ensure rapid heating and a homogeneous and swift sample drying.

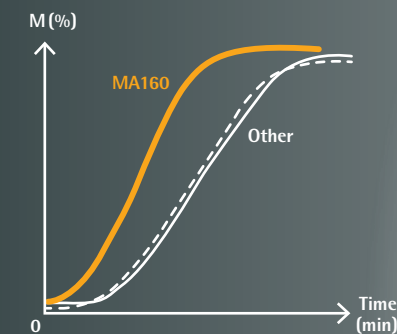


### Large memory capacity

The memory capacity allows the user to store up to 100 different methods and manage them in a method library.

### Fast and precise method development

The MA160's method assistant solves the time-consuming problem of determining and evaluating the right parameters. It enables you to develop new methods quickly in just three simple steps.



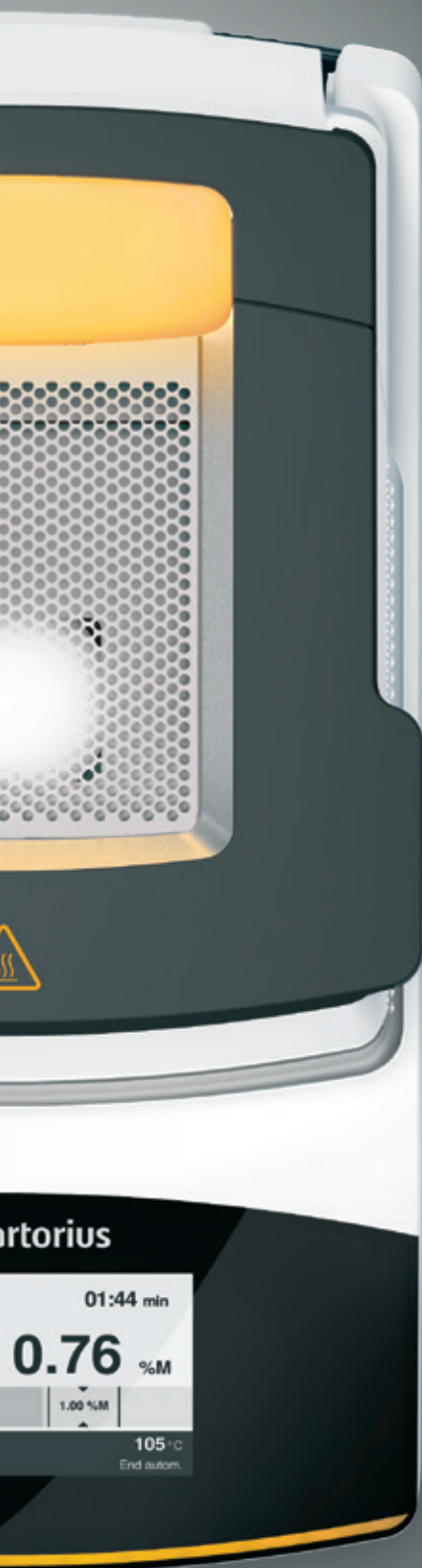

### Display of measurement status

The status lamp indicates the current measurement status during the measurement: "running | START", "process finished | STOP", "analyzer OFF" or "error". It is visible from a distance of up to 10 meters.

### Reliable performance testing

The ReproEasy Pad gives you the option to regularly verify the functionality of the MA160. It guarantees reliable results during routine use.

### User-friendly operation

The intuitive user interface, including touch-screen and easy-to-understand menu, considerably simplifies operation of the MA160.

### Effortless cleaning

BetterClean design enables the user to easily and thoroughly clean the instrument parts. Additionally, the heating module and the sample chamber plate are removable and dishwasher safe.



## Technical Specifications

Max. weighing capacity	200 g
Repeatability, typical	Starting at an initial sample weight of approx. 1 g: $\pm 0.2\%$ , Starting at an initial sample weight of approx. 5 g: $\pm 0.05\%$
Readability	1 mg, 0.01%
Typical sample quantity	5 – 15 g
Reading display	Moisture content in % M and % g   dry matter in % S and g   ATRO in % M/S
Temperature range and settings	40°C–160°C, in increments of 1°C Standby temperature selectable from 40 to 100°C
Sample heating	Infrared heating using an AURI heater, 600 W
Heating programs	Standard drying, gentle drying
Shutoff parameter	Fully automatic, semi-automatic, manual and with timer settings
Sample forceps	Easy-to-handle sample dish
Interface	Mini USB, automatic printer detection, direct transfer to Microsoft® programs
Data transfer	SD card, method import and export function
Housing dimensions (W × D × H)	215 × 400 × 210 mm
Weight	Approx. 6.2 kg

### Accessories

6965542	Disposable sample pans, 80 pcs., aluminum, Ø 90 mm
6906940	Glass fiber pad for analysis of pasty and fatty samples, hard quality, 80 pcs., Ø 90 mm
6906941	Glass fiber pad for analysis of liquid and fatty samples, soft quality, 200 pcs., Ø 90 mm
YHP01MA	ReproEasy pads, 10 pcs. for performance testing to verify the repeatability of the analyzer
YCW512-AC-02	External calibration weight, 100 g (E2) with DKD certificate
YDP40	Standard printer
YDP30	Premium GLP laboratory printer
YCC03-D09	Adapter cable for connecting the YDP20-OCE printer



Disposable sample pans



ReproEasy Pad

# Sales and Service Contacts

For further contacts, visit [www.sartorius.com](http://www.sartorius.com)

## Europe

### Germany

Sartorius Lab Instruments  
GmbH & Co. KG  
Weender Landstrasse 94–108  
37075 Goettingen  
Phone +49.551.308.0  
Fax +49.551.308.3289

### France & Suisse Romande

Sartorius France  
2, rue Antoine Laurent de Lavoisier  
ZA de la Gaudrée  
91410 Dourdan  
Phone +33.1.70.62.50.00  
Fax +33.1.64.59.76.39

### Austria

Sartorius Austria GmbH  
Franzosengraben 12  
1030 Vienna  
Phone +43.1.7965760.0  
Fax +43.1.7965760.24

### Belgium

Sartorius Belgium N.V.  
Leuvensesteenweg, 248/B  
1800 Vilvoorde  
Phone +32.2.756.06.71  
Fax +32.2.253.45.95

### Finland & Baltics

Sartorius Biohit Liquid Handling Oy  
Laippatie 1  
00880 Helsinki  
Phone +358.9.755.951  
Fax +358.9.755.95.200

### Hungary

Sartorius Hungária Kft.  
Kagyló u. 5.  
2092 Budakeszi  
Phone +3623.457.227  
Fax +3623.457.147

### Ireland

Sartorius Ireland Ltd.  
Unit 41, The Business Centre  
Stadium Business Park  
Ballycoolin Road  
Dublin 11  
Phone +353.1.8089050  
Fax +353.1.8089388

### Italy

Sartorius Italy S.r.l.  
Viale A. Casati, 4  
20835 Muggiò (MB)  
Phone +39.039.4659.1  
Fax +39.039.4659.88

### Netherlands

Sartorius Netherlands B.V.  
Phone +31.30.60.53.001  
Fax +31.30.60.52.917  
[info.netherlands@sartorius.com](mailto:info.netherlands@sartorius.com)

### Poland

Sartorius Poland sp.z o.o.  
ul. Wrzesinska 70  
62-025 Kostrzyn  
Phone +48.61.6473830  
Fax +48.61.6473839

### Russian Federation & Ukraine

LLC "Sartorius ICR"  
and LLC "Biohit"  
Uralskaya str. 4, Lit. B  
199155, Saint-Petersburg  
Phone +7.812.327.5.327  
Fax +7.812.327.5.323

### LLC "Biohit"

Post Box 440 "B"  
01001 Kiev, Ukraine  
Phone +380.44.411.4918  
Fax +380.50.623.3162

### Spain & Portugal

Sartorius Spain, S.A.  
Avda. de la Industria, 32  
Edificio PAYMA  
28108 Alcobendas (Madrid)  
Phone Spain +34.902.123.367  
Phone Portugal +351.800.855.800  
Fax Spain +34.91.358.96.23  
Fax Portugal +351.800.855.799

### Switzerland

Sartorius Mechatronics Switzerland AG  
Ringstrasse 24a  
8317 Tagelswangen (ZH)  
Phone +41.44.746.50.00  
Fax +41.44.746.50.50

### U.K.

Sartorius UK Ltd.  
Longmead Business Centre  
Blenheim Road, Epsom  
Surrey KT19 9QQ  
Phone +44.1372.737159  
Fax +44.1372.726171

## America

### USA

Sartorius Corporation  
5 Orville Drive, Suite 200  
Bohemia, NY 11716  
Phone +1.631.254.4249  
Toll-free +1.800.635.2906  
Fax +1.631.254.4253

### Argentina

Sartorius Argentina S.A.  
Int. A. Avalos 4251  
B1605ECS Munro  
Buenos Aires  
Phone +54.11.4721.0505  
Fax +54.11.4762.2333

### Brazil

Sartorius do Brasil Ltda  
Avenida Senador Vergueiro 2962  
São Bernardo do Campo  
CEP 09600-000 - SP- Brasil  
Phone +55.11.4362.8900  
Fax + 55.11.4362.8901

### Canada

Sartorius Canada Inc.  
2179 Dunwin Drive #4  
Mississauga, ON L5L 1X2  
Phone +1.905.569.7977  
Toll-Free +1.800.668.4234  
Fax +1.905.569.7021

### Mexico

Sartorius de México S.A. de C.V.  
Circuito Circunvalación Poniente  
No. 149  
Ciudad Satélite  
53100, Estado de México  
México  
Phone +52.5555.62.1102  
Fax +52.5555.62.2942

## Asia | Pacific

### Australia

Sartorius Australia Pty. Ltd.  
Unit 5, 7-11 Rodeo Drive  
Dandenong South Vic 3175  
Phone +61.3.8762.1800  
Fax +61.3.8762.1828

### China

Sartorius Scientific  
Instruments (Beijing) Co., Ltd.  
33 Yu An Road, Airport Industrial Park  
Zone B,  
Shunyi District, Beijing 101300,  
P.R.China  
Phone +86.10.8042.6300  
Fax +86.10.8042.6486

### Hong Kong

Sartorius Hong Kong Ltd.  
Unit 1012, Lu Plaza  
2 Wing Yip Street  
Kwung Tong  
Kowloon, Hong Kong  
Phone +852.2774.2678  
Fax +852.2766.3526

### India

Sartorius Weighing India Pvt. Ltd.  
#69/2-69/3, NH 48, Jakkasandra,  
Nelamangala Tq  
562 123 Bangalore, India  
Phone +91.80.4350.5250  
Fax +91.80.4350.5253

### Japan

Sartorius Stedim Japan K.K.  
4th Fl., Daiwa Shinagawa North Bldg.  
8-11, Kita-Shinagawa 1-chome  
Shinagawa-ku, Tokyo, 140-0001 Japan  
Phone +81.3.4331.4300  
Fax +81.3.4331.4301

### Malaysia

Sartorius Malaysia Sdn. Bhd  
Lot L3-E-3B, Enterprise 4  
Technology Park Malaysia  
Bukit Jalil  
57000 Kuala Lumpur, Malaysia  
Phone +60.3.8996.0622  
Fax +60.3.8996.0755

### Singapore

Sartorius Singapore Pte. Ltd  
1 Science Park Road,  
The Capricorn, #05-08A,  
Singapore Science Park II  
Singapore 117528  
Phone +65.6872.3966  
Fax +65.6778.2494

### South Korea

Sartorius Korea Ltd.  
8th Floor, Solid Space B/D,  
PanGyoYeok-Ro 220, Bundang-Gu  
SeongNam-Si, GyeongGi-Do, 463-400  
Phone +82.31.622.5700  
Fax +82.31.622.5799

### Thailand

Sartorius (Thailand) Co. Ltd.  
129 Rama 9 Road,  
Huaykwang  
Bangkok 10310  
Phone +66.2643.8361-6  
Fax +66.2643.8367



◀ [www.sartorius.com](http://www.sartorius.com)

Microsoft® is a registered trademark of Microsoft Corporation, USA.